

Preliminary

Product Information Sheet

Microanalytical Reference Material

OKUM*-NP

Nano-particulate pressed powder pellet

Reference Values

Analyte	Value	Uncertainty (95% CL)	Unit	Analyte	Value	Uncertainty (95% CL)	Unit
Na2O	1.136	0.022	g/100g	Sr	16.1	1.0	µg/g
MgO	21.29	0.10	g/100g	Υ	9.08	0.29	µg/g
Al2O3	7.97	0.04	g/100g	Zr	17	1.4	µg/g
SiO2	44.14	0.14	g/100g	Nb	0.37	0.060	µg/g
P2O5	0.0266	0.0023	g/100g	Cs	0.184	0.003	µg/g
K20	0.044	0.002	g/100g	La	0.412	0.017	µg/g
CaO	7.85	0.06	g/100g	Ce	1.27	0.03	μg/g
TiO2	0.38	0.004	g/100g	Pr	0.235	0.008	μg/g
MnO	0.1813	0.0027	g/100g	Nd	1.494	0.06	µg/g
Fe2O3(t)	11.81	0.05	g/100g	Sm	0.715	0.011	µg/g
Sc	27.9	1.5	μg/g	Eu	0.3	0.007	µg/g
V	167.8	3.1	μg/g	Dy	1.61	0.04	µg/g
Cr	2460	31	μg/g	Но	0.355	0.009	µg/g
Со	88.9	1.5	μg/g	Er	1.041	0.014	µg/g
Ni	886	10	μg/g	Yb	1.009	0.023	µg/g
Cu	43.5	1.2	µg/g	Lu	0.148	0.005	μg/g
Zn	61.2	1.9	µg/g	Hf	0.551	0.023	µg/g
Ga	8.79	0.16	µg/g	Ta	0.0264	0.0038	µg/g
₹b	0.96	0.06	µg/g	Th	0.031	0.004	μg/g

Information Values

Analyte	Value	Uncertainty (95% CL)	Unit	Analyte	Value	Uncertainty (95% CL)	Unit
Li	4.4	0.3	μg/g	Gd	1.17	0.07	µg/g
Be	0.065	0.004	μg/g	Tb	0.229	0.010	µg/g
Sn	0.25	0.02	μg/g	Tm	0.155	0.006	µg/g
Sb	0.079	0.012	μg/g	Pb	0.26	0.02	μg/g
Ва	6.2	0.5	µg/g	U	0.012	0.005	μg/g

All values including their uncertainties are taken from the certificate of analysis for the original powder. Information values did not fulfil all necessary statistical criteria of a reference value and should neither be considered for calibration nor validation.

^{*}The original manufacturer (IAG) is not liable for any issues occurring from the use of this material since they took no part in the manufacturing of the pellets.



Intended Use

This microanalytical reference material (MRM) is designed for use by laboratories undertaking the determination of major and trace element mass fractions in komatiite and equivalent matrices with LA-ICP-MS (Laser Ablation Inductively Coupled Plasma Mass Spectrometry), µXRF/XRF (Micro X-ray Fluorescence Spectroscopy) and LIBS (Laser-Induced Breakdown Spectroscopy). It is suitable for calibration and as a secondary reference material for the assessment of a measurement procedure and quality control. Note that the material may only be used for a single purpose in the same measurement process. For example, it must not be used for calibration and method validation at the same time.

Description of the MRM

This MRM is a nanoparticular pressed powder pellet of the komatiite powder "OKUM". The original powder, purchased from the International Association of Geoanalysts (IAG), was subjected to our own material-specific milling protocol and pressed without any binders using a programmable hydraulic press. The fortification of contrasting colour surrounding the reference material is, according to the manufacturer, an "organic compound". The exact composition is not specified any closer. The certificate of analysis is available on demand.

Handling advice and Storage

Avoid touching the pellet's surface directly in order to prevent contamination. Also, do not clean the surface with any liquids as it may compromise the pellet's integrity.

Please note the label marks the bottom of the pellet.

If using a pressed pellet not ordered specifically for μ XRF and or XRF please consider the sample thickness. Store the MRM in a desiccator and or in a dark and dry environment.

The myStandards GmbH cannot be held responsible for changes that happen during storage of the material at the customer's premises, especially with respect to opened samples.

Period of Validity

Provided the storage and handling conditions are met, no chemical alteration is known to exist, and the assigned values will remain stable. Therefore, the product information and assigned values for this MRM are valid for one year from the date of dispatch. This validity may be extended as further evidence of stability becomes available. The manufacturer will inform the customer if any alterations occur.

Date of dispatch: {LIEFERDATUM}

Safety instructions

Nano-particulate powders can cause harm if ingested, inhaled or in contact with skin. In their pressed form however, they do not exhibit any dusting. If a pellet should accidently break, we advise wearing a dust mask during clean up.



Further Information

This MRM has been produced in accordance with the recommendations specified in ISO Guides 30 to 35. Due to processing a part of the sample material may be seen on the fortification, this does not reduce the performance of the MRM. Please refrain from using this part of the pellet.

The pellets are sold exclusively via the myStandards GmbH and authorised subcontractors.

Pellet serial number: {SERIENNUMMER}

Manufactured for: {METHODE}
Size: {GROESSE}

Document History

Version	Date	Changes applied		
1.0	21.01.2019	First publication		
1.1	24.04.2019	24.04.2019 Changed Description of Assigned Values, Description of the MRM, Handling advice and Storage, Period of Validity and Further Information; Insert footnote to Legal notice; Added Document History		
2.0	06.11.2020	Changed Description of the MRM, Handling advice and Storage, Further Information, Intended Use (adding LIBS, XRF and calibration) and Period of Validity (including changing date of manufacturing to date of dispatch); Addition of new terms and conditions; Updated document structure, Addition of individual Pellet characteristics		
3.0	24.06.2022	New official letterhead		
4.0	24.04.2023	Updated link to terms and conditions; Adaptation to automatically in the date and individual pellet characteristics		

References

Potts P.J., Meisel T. Certificate of Analysis, IAG OKUM (Ultramafic rock), International Association of Geoanalysts, 2015, available online at www.iageo.com.

Legal notice

Our order, sales and delivery conditions apply. The valid version of our general terms and conditions (status 01.09.2019) - can be found on our website: https://www.my-standards.com/terms-and-conditions/. They are also available on request.